

PRODUCT INFORMATION

**6X DNA**

**Loading Dye &  
SDS Solution**

#R1151      5 x 1 mL

Lot \_      Expiry Date: \_\_\_\_\_

Store at room temperature or at 4°C for periods up to 12 months. For longer periods, store at -20°C.



**Description**

The 6X DNA Loading Dye & SDS Solution is recommended for preparation of DNA samples with high amounts of DNA binding proteins prior to loading on agarose and polyacrylamide gels. It is supplemented with 1% SDS to eliminate DNA-protein interactions. Proteins have to be denatured to release DNA from DNA-protein complexes. Otherwise, the DNA binding proteins may alter electrophoretic DNA migration, cause DNA samples to stick to gel wells or band shifts. SDS also prevents the annealing of DNA molecules via long cohesive ends and helps to avoid additional bands in electrophoresis. The optimized 6X DNA Loading Dye & SDS Solution contains two electrophoresis tracing dyes: bromophenol blue and xylene cyanol FF. It also contains 100 mM EDTA, that binds divalent metal ions, inhibits metal dependent enzymatic reactions and protects DNA from metal dependent nucleases, effectively inactivating enzymes and protecting DNA from nucleases.



## Composition

0.03% bromophenol blue

0.03% xylene cyanol FF

60% glycerol

1% SDS

100 mM EDTA (pH 7.6, adjusted with Tris)

## Usage Recommendations

- ➊ Add 1 volume of 6X DNA Loading Dye & SDS Solution to 5 volumes of DNA sample.
- ➋ Mix well.
- ➌ Heat at 65°C for 10 minutes.
- ➍ Chill on ice, spin down and load.

## Note

- In 1% agarose gels bromophenol blue co-migrates with ~300 bp DNA, while xylene cyanol FF co-migrates with ~4000 bp DNA.
- The prepared sample can be stored at -20°C. It can be subjected and is viable for several freeze-thaw cycles.

## CERTIFICATE OF ANALYSIS

Tested for DNA sample preparation prior to agarose gel electrophoresis.

Quality authorized by:

 Jurgita Zilinskiene

### LIMITED USE LABEL LICENSE: Internal Research and Development Use Only.

The purchase of this product conveys to the buyer the limited, non-exclusive, non-transferable right (without the right to resell, repackage, or further sublicense) to use this product for internal research and development purposes. No other license is granted to the buyer whether expressly, by implication, by estoppel or otherwise. In particular, the purchase of the product does not include or carry any right or license to use, develop, or otherwise exploit this product commercially and no rights are conveyed to the buyer to use the product or components of the product for purposes including but not limited to provision of services to a third party, generation of commercial databases or clinical diagnostics. This product is sold pursuant to authorization from Thermo Fisher Scientific and Thermo Fisher Scientific reserves all other rights. For information on purchasing a license for uses other than internal research and development purposes, please contact [outlicensing@lifetech.com](mailto:outlicensing@lifetech.com) or Out Licensing, Life Technologies Inc., 5791 Van Allen Way, Carlsbad, California 92008.

### PRODUCT USE LIMITATION

This product is developed, designed and sold exclusively *for research purposes and in vitro use only*. The product was not tested for use in diagnostics or for drug development, nor is it suitable for administration to humans or animals. Please refer to [www.thermoscientific.com/onebio](http://www.thermoscientific.com/onebio) for Material Safety Data Sheet of the product.

© 2015 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries.